

Instructions for Dino-Lite AM / AD 412, 4012 series

Connecting to a television:

It is recommended using an LCD television or Plasma television for better image results. Older generation televisions with video input features will work also. For televisions without direct video input jack, an external device such as a VCR or DVD player with video input can be utilized.

Follow this procedure to connect model AM / AD 412, 4012 series to a television.

IMPORTANT: *The television must have an available video input jack to complete this procedure.*

1. Plug in the AC power adapter into a power outlet (Figure 1).
2. Connect the female power cable end to the supplied male AC power adapter (Figure 2).
3. Connect the composite video cable (yellow) to an available **VIDEO** input jack (yellow) on the television (Figure 3)
4. Turn on the television.
5. Tune the television to the appropriate video input channel to receive the signal from the digital microscope.

NOTE: *If necessary, consult the instruction manual of your television for information about tuning the TV to the correct video input channel.*

How to use:

Simply turn the adjustment dial left or right to adjust focus and magnification.

Included in package

- . Dino-Lite microscope AM / AD 412, 4012 series
- . AC power adapter (110V)
- . Cradle stand

